

RNZO

RNase Decontamination Solution

Cat NO: RN983048, RN983018

Description

RNZO reagent is used to remove RNase contamination from plastic and glass surfaces. It effectively removes high levels of RNase contamination.

Specification

- Store at room temperature (15-25°C).
- Due to chemical material usage that may be hazardous, always make sure to wear suitable lab coat, disposable gloves, and protective eyewear. Material Safety Data Sheet (MSDS) for all products and reagents are provided. They are accessible online at www.rojetechnologies.co.

Before use

Shake or heat at 37°C to bring the precipitate back to solution.

Cleaning work Surface

Apply the RNZO directly to surface to be cleaned, wipe thoroughly with paper towel, rinse with water and then dry with clean paper towel.